




INSTRUCTIONS

Fast Release Pipette Pump™ II

Catalog Number F37911-1002, 1010, 1025

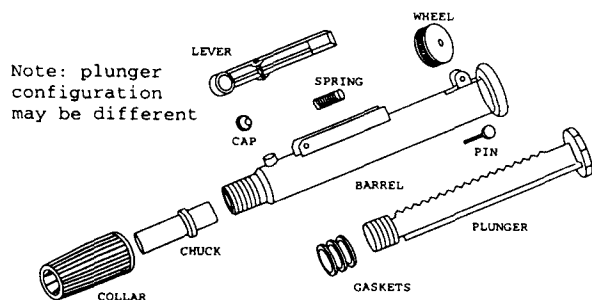

Note: plunger configuration may be different

NOTE: The first time that a pipette is inserted into the pipette pump, the stabilizing fingers at the end of the collar will appear stiff. This is normal for first time use.

Replacement Parts for:

Fast Release Pipette Pump 37911-0000

- 93791-1100 Button
- 93791-1101 Collar
- 93791-1102 Lever
- 93791-1103 Chuck
- 93791-0006 Spring Compression
- 93791-1127 Button & Level Assembly (cap)

Fast Release Pipette Pump 37911-1002 Blue, 2ml

- 93791-1104 Blue Body
- 93789-7001 Rack & Pinion Wheel & Plunger
- 93789-0071 Washer Set

Fast Release Pipette Pump 37911-1010 Green, 10ml

- 93791-1105 Green Body
- 93789-8001 Rack & Pinion Wheel & Plunger
- 93789-0081 Washer Set

Fast Release Pipette Pump 37911-1025 Red, 25ml

- 93791-1106 Red Body
- 93789-9001 Rack & Pinion Wheel & Plunger
- 93789-0091 Washer Set

BEL-ART PRODUCTS®

PEQUANNOCK, NJ 07440-1992
TEL: 973 694-0500 • FAX: 973 694-7199 • www.bel-art.com



INSTRUCTIONS

IMPORTANT NEW PRODUCT INFORMATION

NOTICE OF PRODUCT CHANGE

We are proud to announce the Scienceware® Fast Release Pipette Pump™ II, a new and improved version of our best selling Scienceware® Fast Release Pipette Pump™. The product has been totally redesigned and will replace existing models immediately.

Please use the new catalog numbers listed below when placing future orders for the Scienceware® Fast Release Pipette Pump™ II.

For Pipettes	Replaces Catalog Number	New Catalog Number
Up to 2ml	F37911-0002	F3791-1002
Up to 10ml	F37911-0010	F37911-1010
Up to 25ml	F37911-0025	F37911-1025

The Scienceware® Fast Release Pipette Pump™ II possesses a unique collar that requires no adjustment. The collar clamps the chuck ring to the barrel and holds the chuck in proper alignment, to accept a pipette. Simply push the pipette through the stabilizing fingers at the end of the collar and into the chuck. The stabilizing fingers will prevent the pipette from wobbling while in use.

Insert the pipette into the collar of the pipette pump. Push gently past the stabilizing fingers at the end of the pump's collar. A gentle twisting motion will ease insertion of the pipette. Push the pipette in until it is well seated in the soft chuck.

To pipette-

1. Depress the plunger.
2. Insert proper pipette.
3. Rotate wheel counter clockwise, the liquid will rise in the pipette.
4. Stop when the required amount is reached.
5. To release the liquid drop by drop, turn the knurled wheel slowly in the opposite direction.
6. For fast release of the liquid, press the side lever until the pipette is empty.

WARNING TO ALL GLASS PIPETTE USERS! Caution is advised when using glass pipettes with this product. Glass may break and shatter when forced. Consult your lab safety officer for proper protocol on glass handling.

For disassembly of the pipette pump, please see the diagram at left.

For use with pipettes which retain the last drop.

BEL-ART PRODUCTS®

PEQUANNOCK, NJ 07440-1992
TEL: 973 694-0500 • FAX: 973 694-7199 • www.bel-art.com